
VON MARCKEN
LABORATORY & MEDICAL EQUIPMENT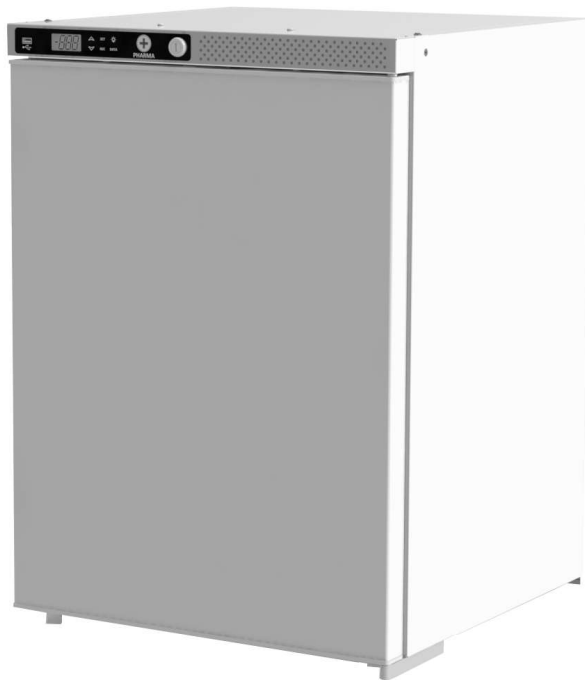

 **AKS 157**

GB Instructions for use.
DE Bedienungsanleitung
FR Mode d'emploi

Warning

As the appliance contains a flammable refrigerant, it is essential to ensure that the refrigerant pipes are not damaged.

The quantity and type of the refrigerant used in your appliance is indicated on the rating plate.

Standard EN378 specifies that the room in which you install your appliance must have a volume of 1m³ per 8 g of hydrocarbon refrigerant used in the appliances. This is to avoid the formation of flammable gas/air mixtures in the room where the appliance is located in the event of a leak in the refrigerant circuit.

WARNING:

Keep ventilation openings in the appliance's cabinet or in the built-in structure clear of obstruction

WARNING:

Do not use other **mechanical devices** or other means to accelerate the defrosting process than those recommended by the manufacturer

WARNING:

Do not damage the refrigerant system

WARNING:

Do not use **electrical appliances** inside the refrigerated storage compartment, unless they are of a type recommended by the manufacturer

WARNING:

Do not expose the appliance to rain, and secure not splashing water when cleaning the floor

WARNING:

This appliance is not intended for use by children or infirm persons unless they have been adequately supervised by a responsible person to ensure that they can use the appliance safely. Children should be supervised to ensure that they do not play with the appliance

WARNING:

If a not compatible USB pen drive is used it can cause a reset of the controller.

**WARNING:**

Do not store explosives, such as aerosol cans with flammable propellants in the unit.

WARNING:

Danger risk of fire or explosion if flammable refrigerant are used. To be repaired only by trained personnel.

WARNING:

Sharp edges on cabinet, compressor compartment, evaporator, ventilation cover and on internal equipment can occur. Please be aware to avoid injury.

WARNING:

The condenser on the back of the appliance will in some cases have a hot surface. Please be aware to avoid injury.

WARNING:

The appliance must be connected to power minimum 12 hours before using it for storage of medicine.

- Always keep the keys in a separate place and out of reach of children
- Do not step on the lower panel to reach medicine in the top of the appliance.
- Before servicing or cleaning the appliance, unplug the appliance from the mains or disconnect the electrical power supply
- If the supply cord is damaged, it must be replaced by the manufacturer, its service agent, or similarly qualified persons in order to avoid a hazard
- **Relevant for Australia:** Supply cord fitted with a plug complies with AS/NZS 3112.
- Frost formation on the interior evaporator wall and upper parts is a natural phenomenon. Therefore, the appliance should be defrosted during normal cleaning or maintenance.
- Please note that changes to the appliance construction will cancel all warranty and product liability


- This device is intended to be used exclusively for medical products.
- If medicine is spilled in the appliance or the defrost water canal is has to be cleaned immediately to avoid the medicine to evaporate to the surroundings.
- If the instructions is lost please contact your supplier of the appliance to have a new instruction for use.
- If service needed to this device, please be aware of only using service personnel with education in handling medical devices.

Contents

Warning	2
Get to know your refrigerator.....	5
Unpacking the appliance.....	6
Before starting the appliance.....	7
Electrial connection.	7
Battery backup system at start-up	8
Installation and start-up	9
Interior fitting.....	10
Maximum shelf and drawer load.....	11
Mounting of drawer dividing system	12
Ventilation guards.....	12
Porthole	13
Reversible door	14
Controller. Operation and function.....	17
Defrosting, cleaning and maintenance	221
Fault finding	22
Warranty, sparepart and service.....	23
Disposal.....	24

Get to know your refrigerator

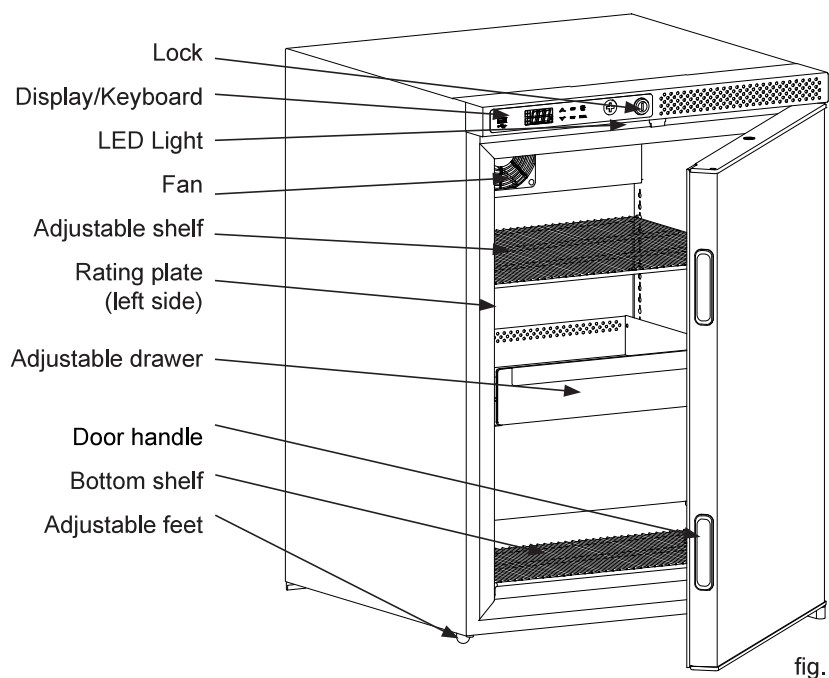Unpacking the appliance

The appliance is delivered in a re-usable wooden box.

1. Reusable wooden package


2. Cut the strapex bands or see next page if sealed with plast clips

3. Remove top

4. Remove top foam inlays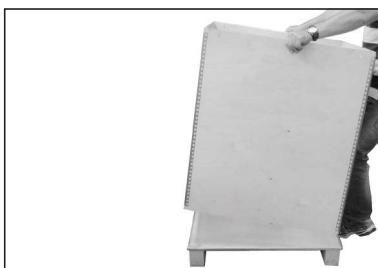

5. Remove box and lift down the appliance

Before starting the appliance

Before operating your new appliance, please read the following instructions carefully as they contain important information on safety, installation, operation and maintenance. Keep the instructions for future reference.

On receipt, check to ensure that the appliance has not been damaged during transport. Transport damage should be reported to the local distributor before the appliance is taken into use.

Clean the inside of the cabinet according to the cleaning instructions.

If during transport the appliance has been laid down, or if it has been stored in cold surroundings (colder than +5°C), it must be allowed to stabilise in an upright position for at least an hour before being switched on.

Note: If the appliance is to be laid down, the door must face upwards and the appliance must be enclosed in the original packaging.

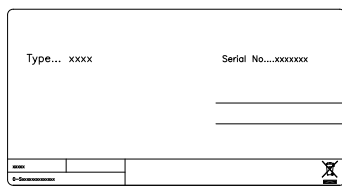
fig. 2

Electrical connection

Wiring and connections in power supply systems must be all applicable (local and national) electrical codes. Consult these codes lengths and sizes prior to cabinet installation.

This device complies with relevant EU directives including Low Voltage Directive 2014/35/EU and Electromagnetic Compatibility Directive 2014/30/EU

The socket should be freely accessible.

Connect the appliance only to 220/240 V / 50Hz alternating current via a correctly installed earthed socket.

The socket must be fused with a 10A or 13A fuse

If the appliance is to be operated in a non-European country, check on the rating plate whether the indicated voltage and current type correspond to the values of your mains supply.

Data regarding voltage and absorbed power / current are given on the rating plate.


The power cord may be replaced by a technician only.

The rating plate provides various technical information as well as type and serial number.


Battery backup system at start-up

This appliance is equipped with a battery back up system.
Please activate the battery backup before final placement (see fig. 3)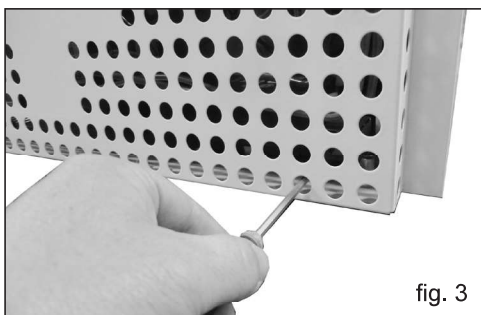

Use a pen or screwdriver to push the button through the above hole on the appliance backside. This will switch on the battery backup system.

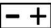Then please check the battery backup image  appears in the front display lower left corner. This shows the battery backup is activated.

NOTE!

The battery backup system does not supply the cooling system with power.

When starting up the appliance for the first time it is necessary to switch on the battery backup system.

NOTE!

The battery for back up should be changed every third year to secure 48 hours of back up. Please put this change in the maintenance schedule for every third year.

Battery backup function

The battery backup system supplies the controller and keyboard with power at power failure. This makes it possible to supervise the temperatures in the unit during the power failure. The battery backup system makes it possible to supervise the temperatures for 48 hours. After a power failure and at the first start up the battery needs to be recharged. To regain the full capacity the battery will be reloading for 10 days

Installation and start-up

The appliance viewed from above

If the appliance is placed beside a wall, there must be sufficient room for its door to be opened wide enough to allow the shelf to be pulled out (fig. 4)

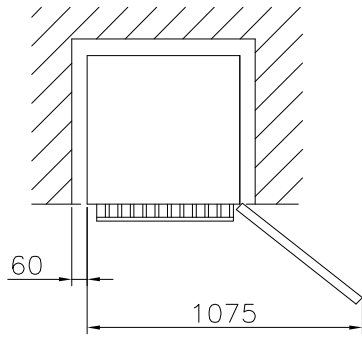
fig. 4

Side by side

In side-by-side arrangements there must be at least 60 mm between each unit on each side of the cabinets, to ensure the doors can be opened freely (See fig. 5)

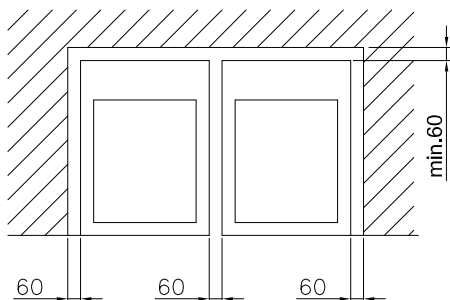
fig. 5

Setting up

It is important that the appliance is absolutely level. It can be levelled by screwing the adjustable feet of the appliance up or down.

Use a spirit level to check that the appliance is absolutely level sideways.

If the appliance is placed on a soft surface, e.g. floorboards or a carpet you must recheck whether the appliance is still level after a period of time as the underlying surface may yield under the weight of the appliance.

Ventilation

It is important that the appliance is well ventilated and that air can circulate freely above, below and around it.

The figures (fig. 4-6) illustrate how the necessary air circulation around the appliance is ensured.

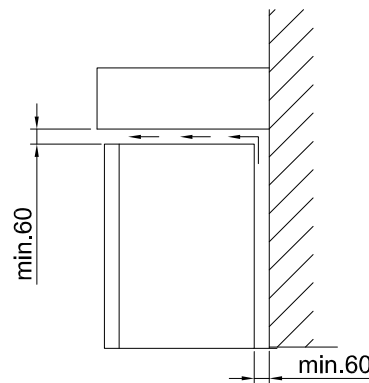
fig. 6

Interior fitting

Bottom wire shelf

The small wire shelf is easily removed & relocated during cleaning or else

Upper wire shelves

The big wire shelves can be changed in height position by following this instruction


1. demount the two fixation clips in the right side
2. slide the shelf, direction right
3. pull up the left side (clips still mounted)
4. push the shelf into the right hole pair for the new wanted shelf height position
5. level the shelf horizontally and pull it into the left hole pair
6. remount the two fixation clips in the right side



Rails & drawers

This instruction tells how to change a drawer height position

1. lift the drawer front and pull outwards until the drawer is free from the rails
2. remove the rails: start by lifting up the closest rail end and pull it out. Then the far rail end is set free by carefully pulling towards yourself
3. remount the rails in the new height position
4. reposition the drawer by carefully lowering it's wheels into both rails. In the same movement push the drawer in final position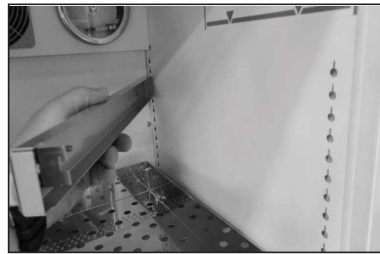


Maximum shelf & drawer load

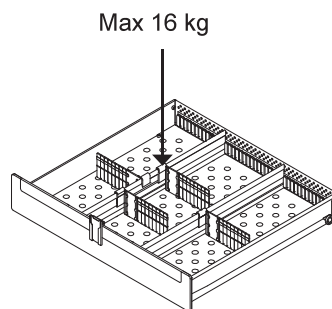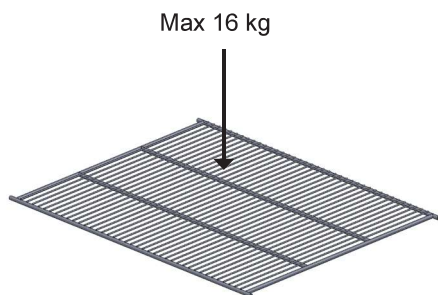
Mounting of drawer dividing system


Drawers are optional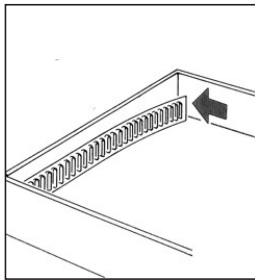

1. The Rail Holder with double adhesive tape is mounted inside the rear and front of the drawer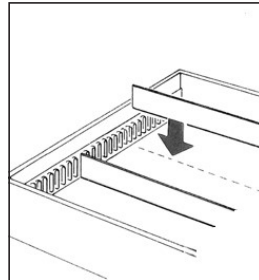

2. Put in the rails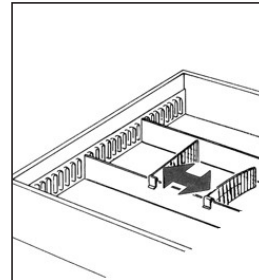

3. Snap on the dividers and slide it

Adjustment of dividing parts

Integrated breakpoints allows storage of individual products. Adjust the dividing part as follows.

(See fig 7) Cut into the appropriate seam and break of the excess part.

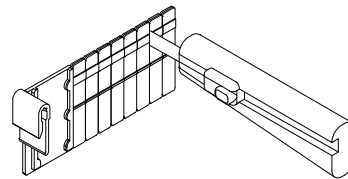
fig. 7

Ventilation guards

The two ventilation guards are mounted before moving the appliance to its final placement. (See fig.8)

The guards function is to secure an absolute minimum of ventilation, if the appliance is pushed too close up against a wall. Remember our guideline is minimum 60 mm free space behind the appliance to ensure sufficient airflow.


The Ø15 mounting holes are located one in each side of the engine cover on the appliance back. The ventilation guards are pushed into the holes


fig. 8

Porthole


The AKG/AKS157 Undercounter has a porthole which can be used for an additional sensor/probe.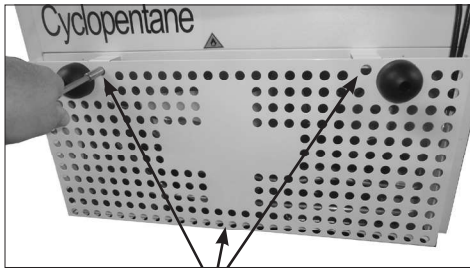


1. Loosen these three screws on the appliance back and remove the cover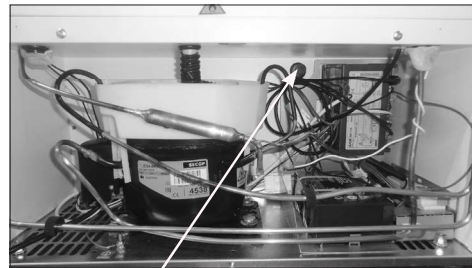

2. The porthole is located on the rear wall (follow arro) Remove the rubber plug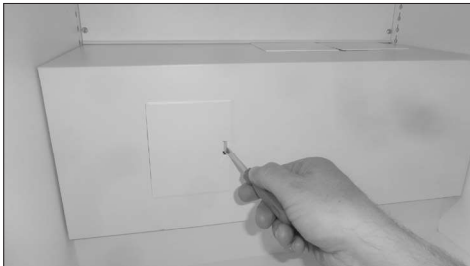

3. Remove this service hatch inside the appliance with a screwdriver

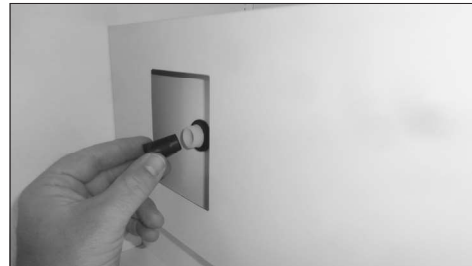4. Remove the rubber plug and guide the sensor/probe through the porthole

5. Re-seal both porthole ends with putty and remount service hatch (inside) + back-cover

GB

Reversible door


1. Switch off main power and place the appliance on a table

2. Open door and remove two screws in top panel


3. Remove lock pin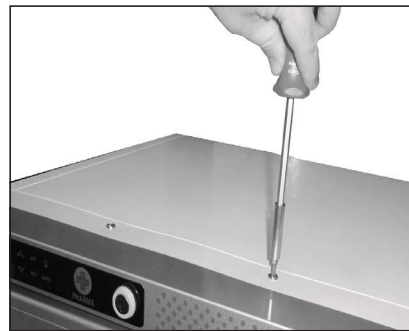


4. Remove four screws from top panel

5. Pull out the top panel and ...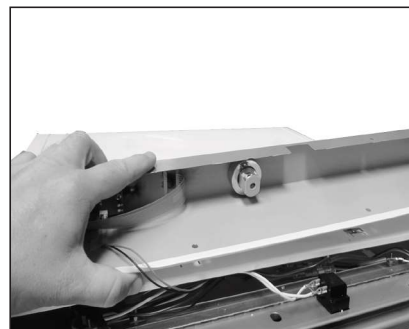


6. ...place it upside down on the cabinet top

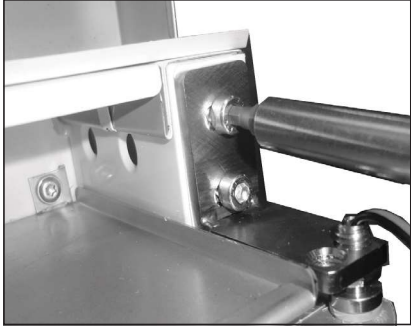7. Remove the two top hinge screws and support the door from tilting afterwards

8. Unscrew the top hinge pivot with a spanner

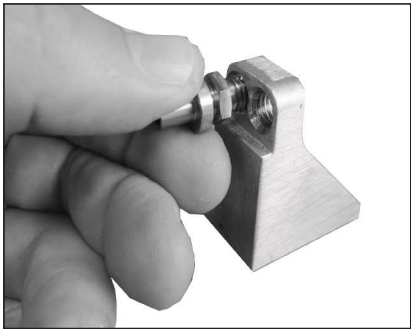9. Mount the pivot in the other hole and tighten it.

10. Gain access below bottom hinge, unscrew foot and...

11. ...demount screws. Then remove bottom hinge and remount in new hinge side when ready


12. In the opposite side: demount the foot + screws...


GB

13. ...and remount them in the former hinge side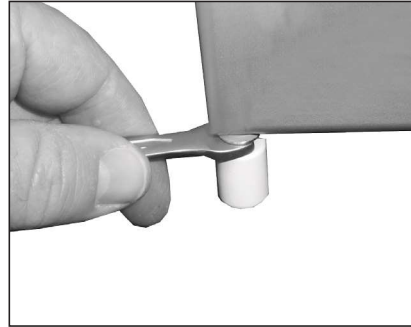

14. Place the door over the bottom hinge while holding the door closer tight. Lower the door pivot carefully down into the hinge

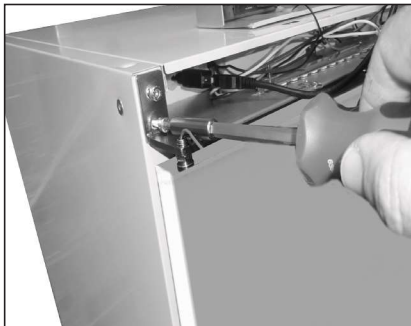15. Remount the top hinge. Make sure the door is centered on the cabinet front before tightening the screws

16. Put the toppanel back in place

17. Remount four screws into toppanel

18. Remount two screws and lock pin in toppanel and the door hinge side change is complete

Controller. Operation and function

User interface

LEDS

Each LED function is described in the following table.

LED	MODE	Function
	ON	Compressor enabled
	Flashing	Anti-short cycle delay enabled
	ON	Defrost enabled
	Flashing	Drip time in progress
	ON	Fans enabled
	Flashing	Fans delay after defrost in progress
	ON	An alarm is occurring
	ON	Recording activated
	ON	Battery is fully charged
	Flashing	Battery is beeing charged
	Flashing	Charging problem or battery failure
°C/°F	ON	Measurement unit
°C/°F	Flashing	Programming phase

Buttons

SET	To display target set point; in programming mode it selects a parameter or confirm an operation.
	To enter fast access menu In programming mode it browses the parameter codes or increases the displayed value. (DOWN); n programming mode
	In programming mode it browses the parameter codes or decreases the displayed value. Push it for 3s to start a manual defrost
DATA	Export data from button
	To switch off the light in the door (only models with light in the door)
REC	Log activation and deactivation from button (Password protected)

How to see and modify the setpoint

How to: See the Set point

1. Push and immediately release the SET key: the display will show the Set point value.
2. Push and immediately release the SET key or wait for 60 sec to display the temperature in the unit

How to: Change the Set point

1. Push the SET key more than 2 sec to change the Set point value.
2. The value of the set point will be displayed and the "°C" LED will start blinking.
3. To change the Set value push the UP or DOWN arrows within 60 sec.
4. To save the new set point value, push the SET key again and wait for 3 sec.

NOTE: To exit without making any change to the set point, push the SET key or wait 60 sec.

Clock settings and RTC alarm reset

1. Push the UP key once, to acces the menu.
2. The display shows H0, then push SET.
3. The parameters for setting time and date occurs. To set the parameter push SET, push the UP or DOWN button to change the parameter. Confirm by pushing SET.

The time and date parameters:

Hur: hour
0in: minutes
Udy: weekday
dAy: date
0on: month
yEA: year

To exit: Press **SET + UP** keys for about 10 sec. or wait 60 sec.

How to export data and alarms to USB

1. Insert the **USB** key
2. Push the **DATA** key for more that 3 sec.
3. Controller starts uploading data to **USB**
4. At the end the following message will be displayed:
 - a. **“End”** if everything is ok
 - b. **“Err”** if exporting has not taken place.

Alarms

The controller memorizes the last 100 alarms happened, together with their start and finish time. It's possible to export the alarms as described in the previous chapter.

Active alarm

Controller will show active alarm alternated with the temperature inside the unit.

Alarm signals

Message	Cause
"PF1"	Regulating probe TR3 failure
"PF2"	EVP Probe Failure
"PF3"	Logging probe TL1 failure
"PF4"	Logging probe TL2 failure
"HA1"	TR3 High Alarm
"LA1"	TR3 Low Alarm
"HA3"	High temperature alarm probe TL1
"LA3"	Low temperature alarm probe TL1
"HA4"	High temperature alarm probe TL2
"LA4"	Low temperature alarm probe TL2
"dA"	Door Open Alarm
"CA"	Serious Alarm

The alarm message is displayed until the alarm condition is recovered.
All the alarm messages are showed alternating with the temperature in the unit.
Except for the "PF1" which is flashing.

Defrosting, cleaning and maintenance

Automatic defrosting.

The refrigerator is defrosted automatically. Defrost water runs through a pipe and is collected in a tray above the compressor where the heat generated by the compressor causes it to evaporate. The defrost water tray should be cleaned when needed or at least once a year

Cleaning and maintenance

Before cleaning, disconnect the power supply and remove all loose fittings.

The appliance is best cleaned using warm water (max. 65°C) with a little mild, perfume-free detergent. Never use cleaning agents that scour. Use a soft cloth. Rinse with clean water and dry thoroughly. It is important to prevent water from entering the control panel.

Add a few drops of disinfectant to the defrost water drain a couple of times a year, and clean the drain using a pipe cleaner or similar. Never use sharp or pointed implements.

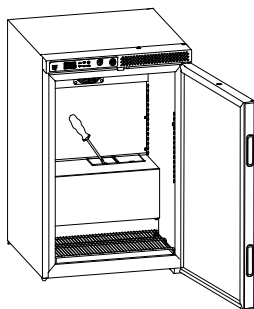
fig.9

Access to the defrost water drain is gained by lifting up the centerplaced servicehatch inside the cooler (See fig. 9)

When disinfecting the appliance always consult professional persons to ensure that the disinfectant used is approved in your country, and to ensure right method is used.

For use in this appliance, we recommend the following base of disinfectant: alcohols and aldehydes. Chlorine and peracids can also be used, but please be aware that chlorine and peracids based products can inflict on the surface of materials, there for a careful washing off with clear water and wiping off, after disinfecting, is very important.

The sealing strip around the door must be cleaned regularly, at least twice a year. to prevent discolouration and prolong service life. Use clean water. After cleaning the sealing strip, check that it continues to provide a tight seal.

It is recommended that the wire and tubular condensers on the rear of the appliance are cleaned.

Dust collecting on the condenser on the rear of the cabinet, the compressor and in the compressor compartment is best removed using a vacuum cleaner. This must be done regularly, at least four times a year

OBS: If the appliance is not to be used for any length of time, switch it off, disconnect the power supply, empty it, clean the inside and leave the door open to allow air circulation and prevent smells.

WARNING: Secure not splashing water when cleaning the floor.

Fault finding

Fault	Possible cause	Remedy
Screen of keyboard is not lit.	Power failure; the fuse is blown; the appliance is not plugged in correctly, the power is cut.	Check that power is connected. Reset the fuse
Temperature in refrigerator too high	The ventilation grille is blocked. The door is not closed properly. The quantity of items placed in the freezer at a time is greater than the capacity of the appliance. The temperature setting is too high	Ensure unhindered air circulation Close the door and wait 15 min. Wait 15min Lower the temperature setting.
Temperature in refrigerator too low.	The temperature setting is too low	Raise the temperature setting
Vibration or bothersome noise	The appliance is not level.	Level the appliance (For model with adjustable feet: use a spirit level)
Compressor runs continuously.	The temperature setting is too low. Too high room temperature.	Raise the temperature setting. Ensure adequate ventilation
Display or alarm log shows "RTC" error	The date and time is not set.	Please set the date and time.

Warranty, spare parts and service

Warranty disclaimer

Faults and damage caused directly or indirectly by incorrect operation, misuse, insufficient maintenance, incorrect building, installation or mains connection. Fire, accident, lightening, voltage variation or other electrical interference, including defective fuses or faults in mains installations.

Repairs performed by others than approved service centres and any other faults and damage that the manufacturer can substantiate are caused by reasons other than manufacturing or material faults are not covered by the warranty.


Please note that changes to the construction of the appliance or changes to the component equipment of the appliance will invalidate warranty and product liability, and the appliance cannot be used lawfully. The approval stated on rating plate will also be invalidated.

Transport damage discovered by the buyer is primarily a matter to be settled between the buyer and the distributor, i.e. the distributor must ensure that such complaints are resolved to the buyer's satisfaction.

Before calling for technical assistance, please check whether you are able to rectify the fault yourself. If your request for assistance is unwarranted, e.g. if the appliance has failed as a result of a blown fuse or incorrect operation, you will be charged the costs incurred by your call for technical assistance.

Spare parts

When ordering spare parts, please state the type, serial and product numbers of your appliance. This information is given on the rating plate. The rating plate contains various technical information, including product-type and serial numbers.

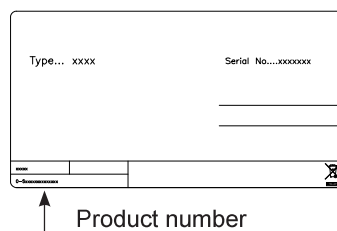Disposal

Information for Users on Collection and Disposal of Old Equipment and used Batteries

These symbols on the products, packaging, and/or accompanying documents mean that used electrical and electronic products and batteries should not be mixed with general household waste. For proper treatment, recovery and recycling of old products and used batteries, please take them to applicable collection points, in accordance with your national legislation and the Directives 2002/96/EC and 2006/66/EC.

By disposing of these products and batteries correctly, you will help to save valuable resources and prevent any potential negative effects on human health and the environment which could otherwise arise from inappropriate waste handling.

For more information about collection and recycling of old products and batteries, please contact your local municipality, your waste disposal service or the point of sale where you purchased the items.

Penalties may be applicable for incorrect disposal of this waste, in accordance with national legislation.

For business users in the European Union.

If you wish to discard electrical and electronic equipment, please contact your dealer or supplier for further information.

Information on Disposal in other Countries outside the European Union

These symbols are only valid in the European Union. If you wish to discard this product, please contact your local authorities or dealer and ask for the correct method of disposal.

Note for the battery symbol:

This symbol might be used in combination with a chemical symbol. In this case it complies with the requirement set by the Directive for the chemical involved.

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A/S Vestfrost